

Registro PCR multisegmento para Influenza A

Zhou B et al. J Virol. 2009 Oct;83(19):10309-13

PrimersUni12/Inf-1 5'- GGG GGG AGC **AAA** AGC AGG -3'Uni12/Inf-3 5'- GGG GGG AGC **GAA** AGC AGG -3'

Uni13/Inf-3 5'- CGG GTT ATT AGT AGA AAC AAG G -3'

Superscript III/Platinum Taq HiFi Enzyme Mix (Invitrogen)

Preparación de la mezcla

Lote	Reactivo	1x (μl)	x (μl)	[] final
[] _____	H2O	10		
[] _____	2X RT-PCR Buffer	25		
[] _____	Uni12/Inf1 10μM	0,8		0,16 μM
[] _____	Uni12/Inf3 10μM	1,2		0,24 μM
[] _____	Uni13/Inf3 10μM	2		0,4 μM
[] _____	SSIII TaqHiFi	1		

PROGRAMA A MULTISE

RT	PCR		
42°C 60'	94°C 30"	94°C 30"	68°C 5'
94°C 2'	44°C 30"	57°C 30"	
	68°C 3'	68°C 3'	
	5x	35x	

Alicuotado y adición del extractoPipetas: P10 ☐ P20 ☐**Termociclador, Base, Modulo Óptico**CC010388 ☐ CT055496 ☐ 785BR 11404 ☐ 03495 ☐ 02472 ☐OPUS 795BR02958 ☐ OPUS 795BR02587 ☐

Añadir 10 μl de extracto a los 40 μl de Mezcla

Revisado por _____

Nº	Resultado	Comentario	Geles
1			
2			
3			
4			
5			
6			
7			
8			
9			
10			
11			
12			
13			
14			
15			
16			
17			
18			
19			
20			